

siRNA Cartridge Purification - Crude Fraction

Column: Clarity® 3 µm Oligo-RP, LC Column 50 x 4.6 mm, Ea

Dimensions: 50 x 4.6 mm ID

Order No: 00B-4441-E0

Elution Type: Gradient

Eluent A: 50mM TEAA, pH 7.5 / 5% ACN

Eluent B: Methanol

Gradient Profile:	Step No.	Time (min)	Pct A	Pct B
	1	0	100	0
	2	20	40	60

Flow Rate: 1 mL/min

Col. Temp.: ambient

Detection: UV-Vis Abs.-Variable Wave.(UV) @ 260 nm (ambient)

Analyst Note: SecurityGuard™ Guard Cartridge System extends column lifetime.


- SecurityGuard Cartridges, Clarity Oligo-RP 4 x 3.0mm, 10/Pk Part No.: AJ0-8135

- Holder Part No.: KJ0-4282

15800



Products used in this application:

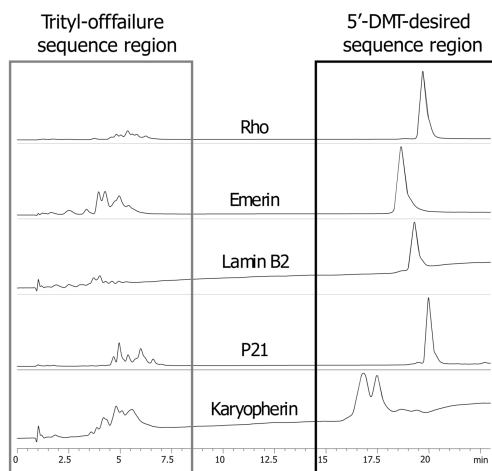

ANALYTES:

1 Crude siRNA

Retention Time: 14 min

