


**Solvents on ZB-WAXplus 30x0.25x0.25****Column:** Zebtron™ ZB-WAXplus™, GC Cap. Column 30 m x 0.25 mm x 0.25 µm, Ea**Phase:** Polyethylene Glycol (PEG)**Dimensions:** 30 meters x 0.25 mm x 0.25 µm**Order No:** 7HG-G013-11**Oven Profile:** 30°C for 1min to 70°C at 14°C/min to 220°C at 25°C/min for 3 min**Carrier Gas:** Constant Flow Helium, 1.2 mL/min**Injection:** Split 50:1 0.2 µL @ 220°C**Detection:** Mass Selective (MSD) (250°C)**Analyst Note:** All analytes are 85 ppm in DMA**Products used in this application:**

**Solvents on ZB-WAXplus 30x0.25x0.25****ANALYTES:**

|                                                     |                                 |
|-----------------------------------------------------|---------------------------------|
| <b>1</b> air                                        | <b>41</b> Butanol               |
| <b>2</b> Pentane                                    | <b>42</b> Nitromethane          |
| <b>3</b> Hexane                                     | <b>43</b> Cumene                |
| <b>4</b> Ethyl ether                                | <b>44</b> 2-Methoxy ethanol     |
| <b>5</b> 2-Methoxy-2-methyl-2-propane               | <b>45</b> Pyridine              |
| <b>6</b> Heptane                                    | <b>46</b> o-Xylene              |
| <b>7</b> Cyclohexane                                | <b>47</b> 3-Methyl butanol      |
| <b>8</b> 1,1-Dichloroethene                         | <b>48</b> Chlorobenzene         |
| <b>9</b> Methylcyclohexane                          | <b>49</b> n-Pentanol            |
| <b>10</b> Acetone                                   | <b>50</b> DMF                   |
| <b>11</b> Formic acid ethyl ester                   | <b>51</b> Anisole               |
| <b>12</b> Acetic acid methyl ester (methyl acetate) | <b>52</b> Tetralin              |
| <b>13</b> Tetrahydrofuran                           | <b>53</b> Dimethyl sulfoxide    |
| <b>14</b> Carbon tetrachloride                      | <b>54</b> Ethylene glycol       |
| <b>15</b> 1,1,1-Trichloroethane                     | <b>55</b> Methylacetamide       |
| <b>16</b> Ethyl acetate                             | <b>56</b> N-Methylpyrrolidinone |
| <b>17</b> Acetic acid-1-methyl ethyl ester          | <b>57</b> Formamide             |
| <b>18</b> 2-Butanone                                |                                 |
| <b>19</b> 1,2-Dimethoxyethane                       |                                 |
| <b>20</b> Dichloromethane                           |                                 |
| <b>21</b> Isopropanol                               |                                 |
| <b>22</b> Benzene                                   |                                 |
| <b>23</b> Propyl acetate                            |                                 |
| <b>24</b> 1,2-Dichloroethene                        |                                 |
| <b>25</b> Trichloroethylene                         |                                 |
| <b>26</b> Acetonitrile                              |                                 |
| <b>27</b> Methylisobutylketone (MIBK)               |                                 |
| <b>28</b> Isobutyl acetate                          |                                 |
| <b>29</b> Chloroform                                |                                 |
| <b>30</b> Toluene                                   |                                 |
| <b>31</b> n-Propanol                                |                                 |
| <b>32</b> Water                                     |                                 |
| <b>33</b> 1,4-Dioxane                               |                                 |
| <b>34</b> 1,2-Dichloroethane                        |                                 |
| <b>35</b> Butyl acetate                             |                                 |
| <b>36</b> 2-Hexanone                                |                                 |
| <b>37</b> 2-Methylpropanol                          |                                 |
| <b>38</b> Ethylbenzene                              |                                 |
| <b>39</b> p-Xylene                                  |                                 |
| <b>40</b> m-Xylene                                  |                                 |

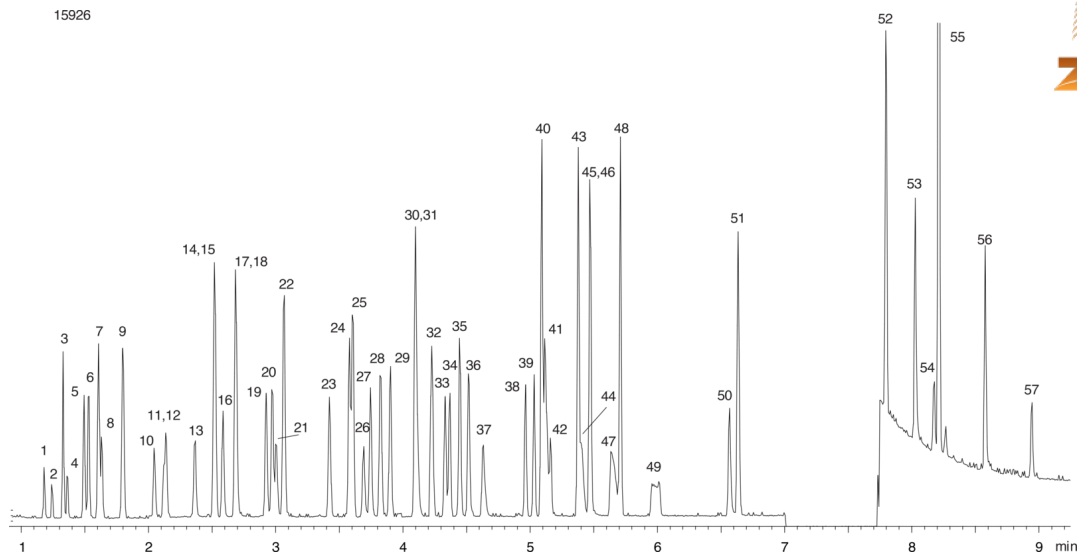