

Amphetamines by MSD on a ZB-5MSi**Column:** Zebtron™ ZB-5MSi, GC Cap. Column 10 m x 0.18 mm x 0.18 µm, Ea**Phase:** 5% Phenyl 95% Dimethylpolysiloxane**Dimensions:** 10 meters x 0.18 mm x 0.18 µm**Order No:** 7CD-G018-08**Oven Profile:** 100 °C to 320 °C @ 25 °C/min**Carrier Gas:** Constant Flow Helium, 1 mL/min**Injection:** Split 10:1 1 µL @ 280°C**Detection:** Mass Selective (MSD) (320°C)**Analyst Note:** Analytes are 40ppm derivatized with HFBA

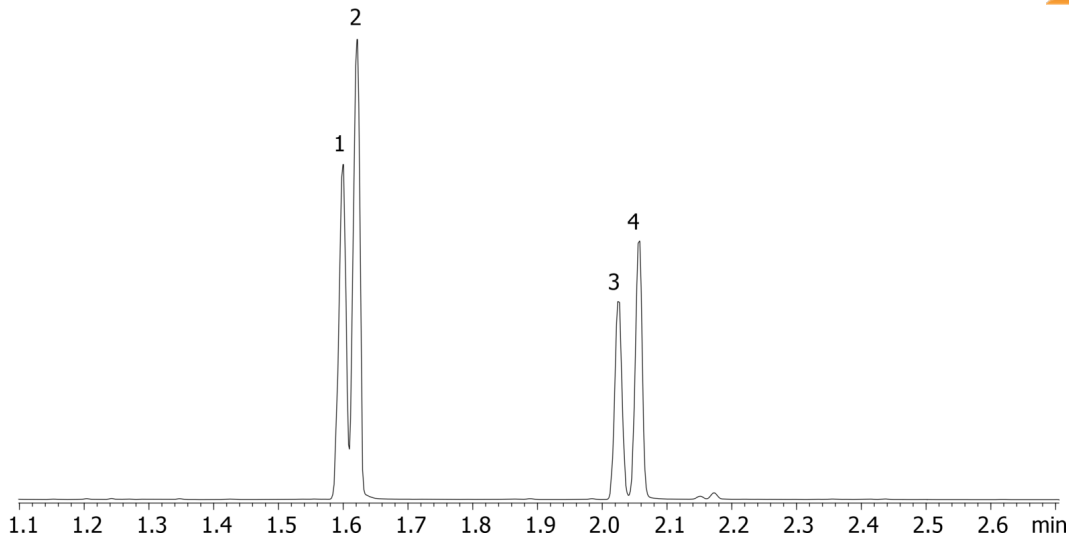
Sim ions:

(Quant ion*) D11-Amphetamine (*244,128): Amphetamine (*240,118,117): D14-Methamphetamine (*261,213):

Methamphetamine (*254,210,118)

**Products used in this application:**

17793

**ANALYTES:**

- 1 Amphetamine-d11
- 2 Amphetamine
- 3 Methamphetamine-d14
- 4 Methamphetamine

