


Underivatized Active Drugs on ZB-XLB-HT**Column:** Zebron™ ZB-XLB-HT, GC Cap. Column 30 m x 0.25 mm x 0.25 µm, Ea**Phase:** Proprietary XLB Phase**Dimensions:** 30 meters x 0.25 mm x 0.25 µm**Order No:** 7HG-G024-11**Oven Profile:** 135 °C to 220 °C @ 10 °C/min to 320 @ 25 °C/min for 8 min**Carrier Gas:** Constant Flow Hydrogen, 40 cm/sec**Injection:** Split 42:1 2 µL @ 250°C**Detection:** Flame Ionization (FID) (325°C)**Analyst Note:** All analytes are 50ppm in methanol

17896

**Products used in this application:****ANALYTES:**

- 1 Ibuprofen
- 2 Dimenhydrinate
- 3 Glutethimide
- 4 Desipramine
- 5 Oxymorphone
- 6 Quinidine
- 7 Alprazolam

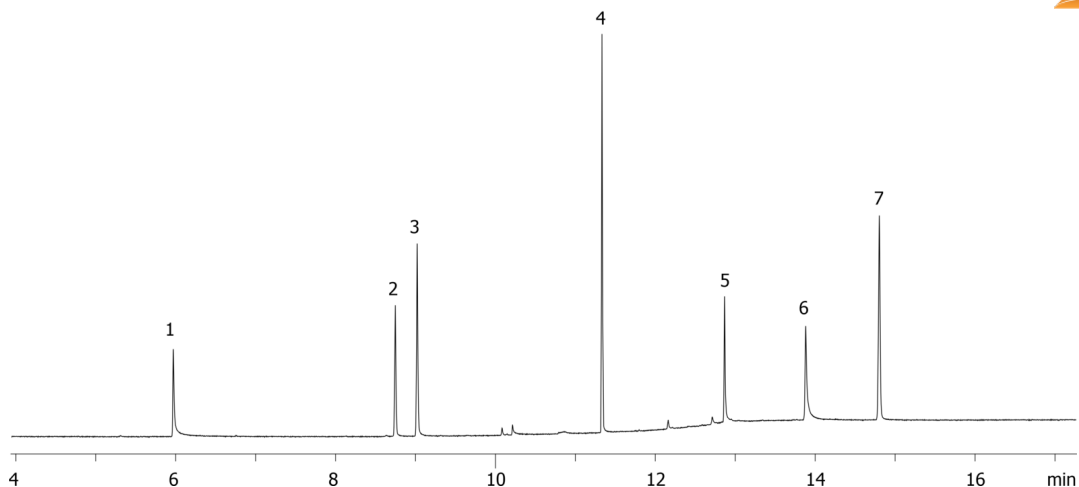