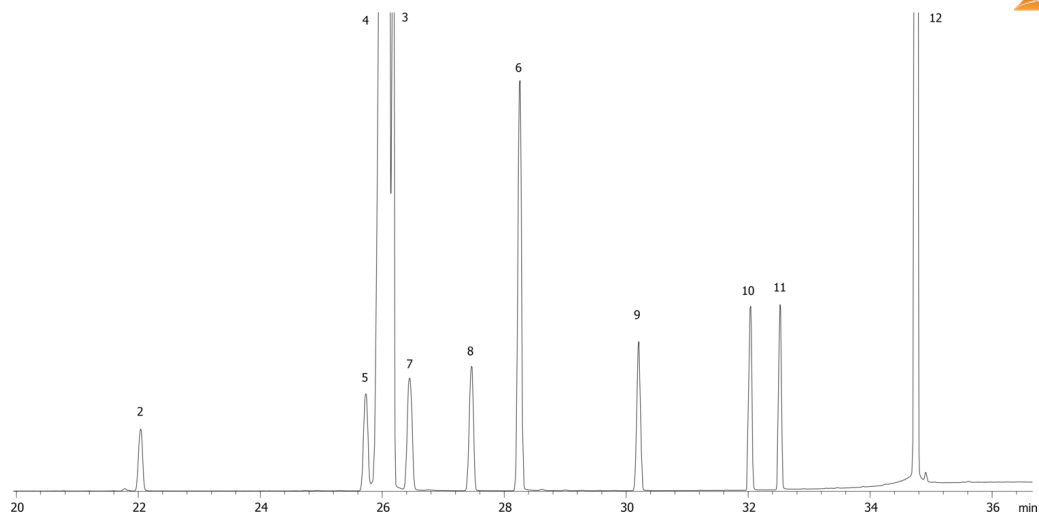


Organohalide Pesticides (EPA 505) Mix # 1 on ZB-1701P

Column: Zebtron™ ZB-1701, GC Cap. Column 30 m x 0.25 mm x 0.25 µm, Ea
Phase: 14% Cyanopropylphenyl 86% Dimethylpolysiloxane
Dimensions: 30 meters x 0.25 mm x 0.25 µm
Order No: 7HG-G006-11
Oven Profile: 40 °C for 1 min to 100 °C @ 15 °C/min to 225 °C @ 5 °C/min to 290 °C @ 15 °C/min for 25 min
Carrier Gas: Constant Flow Helium, 0.8 mL/min
Injection: On-Column :1 0.5 µL @ 43°C
Detection: Mass Selective (MSD) (335°C)
Analyst Note: Analytes 10 ppm in Acetone

**Products used in this application:**

Organohalide Pesticides (EPA 505) Mix # 1 on ZB-1701P

ANALYTES:

- 1 Hexachlorocyclopentadiene
- 2 Hexachlorobenzene
- 3 Simazine
- 4 Atrazine
- 5 gama-BHC
- 6 Alachlor
- 7 Heptachlor
- 8 Aldrin
- 9 Heptachlor epoxide
- 10 Dieldrin
- 11 Endrin
- 12 Methoxychlor

