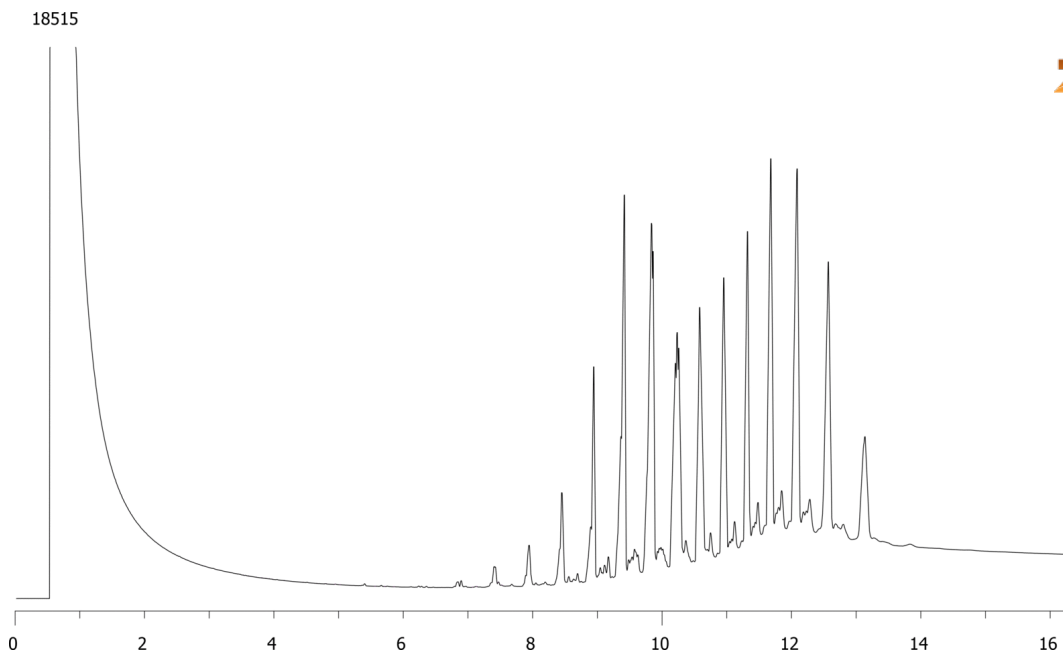


Butter Triglycerides on ZB-XLB-HT by FID**Column:** Zebtron™ ZB-XLB-HT, GC Cap. Column 15 m x 0.25 mm x 0.10 µm, Ea**Phase:** Proprietary XLB Phase**Dimensions:** 15 meters x 0.25 mm x 0.1 µm**Order No:** 7EG-G024-02**Oven Profile:** 150 °C to 380 °C @ 20 °C/min for 5 min**Carrier Gas:** Constant Flow Helium, 1.3 mL/min**Injection:** On-Column :1 1 µL @ 153°C**Detection:** Flame Ionization (FID) (350°C)**Analyst Note:** Butter was 1.0mg/mL of isooctane**Products used in this application:****ANALYTES:****1** butter