

Hu IgG2k aggregates on BioSep3000(2)

Column: BioSep™ 5 µm SEC-s3000 290 Å, LC Column 300 x 7.8 mm, Ea

Dimensions: 300 x 7.8 mm ID

Order No: 00H-2146-K0

Elution Type: Isocratic

Eluent A: 50mM Sodium Phosphate, 300mM NaCL pH 6.8

Gradient	Step No.	Time (min)	Pct A
Profile:	1	0	100

Flow Rate: 1 mL/min


Col. Temp.: ambient

Detection: UV-Vis Abs.-Variable Wave.(UV) @ 220 nm (ambient)


Analyst Note: SecurityGuard™ Guard Cartridge System extends column lifetime.

- SecurityGuard Cartridges, GFC 3000 4 x 3.0mm ID, 10/Pk Part No.: AJ0-4488

- Holder Part No.: KJ0-4282

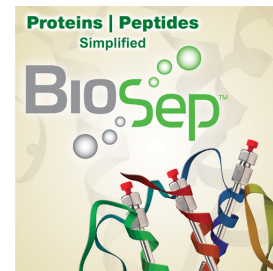

Products used in this application:

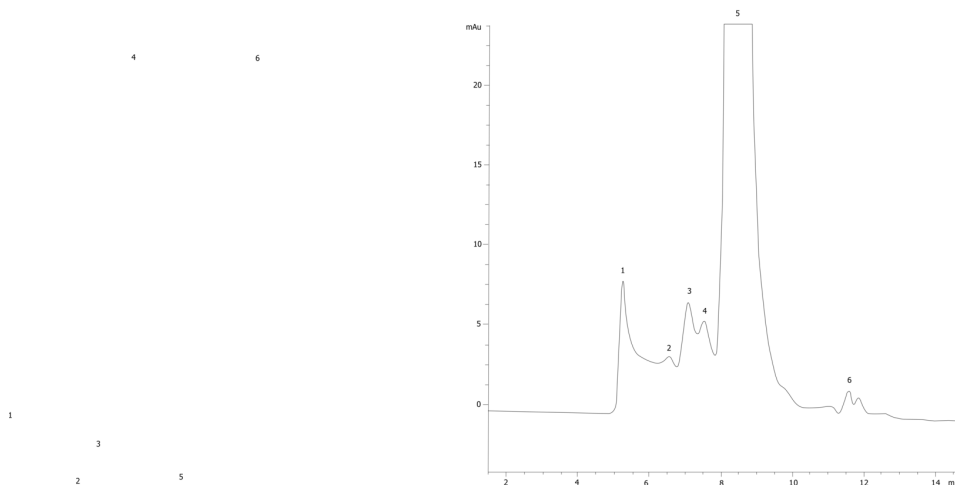

ANALYTES:

- 1 High MW aggregate
Retention Time: 5.26 min
- 2 IgG trimer
- 3 IgG2 dimer
- 4 IgG dimer
- 5 Human IgGk monomer
- 6 void peak

