

Overlay: Yarra 3µm SEC-3000 vs TSKgel G3000 SWxl Low Salt Buffer Competitive Comparisons

Column: Yarra™ 3 µm SEC-3000, LC Column 300 x 7.8 mm, Ea

Dimensions: 300 x 7.8 mm ID

Order No: 00H-4513-K0

Elution Type: Isocratic


Eluent A: 100 mM Sodium Phosphate in Water pH 6.8 with 0.025% Na₃

Gradient Profile:	Step No.	Time (min)	Pct A
1	1	0	100

Flow Rate: 1 mL/min

Col. Temp.: ambient

Detection: UV-Vis Abs.-Variable Wave.(UV) @ 280 nm (ambient)

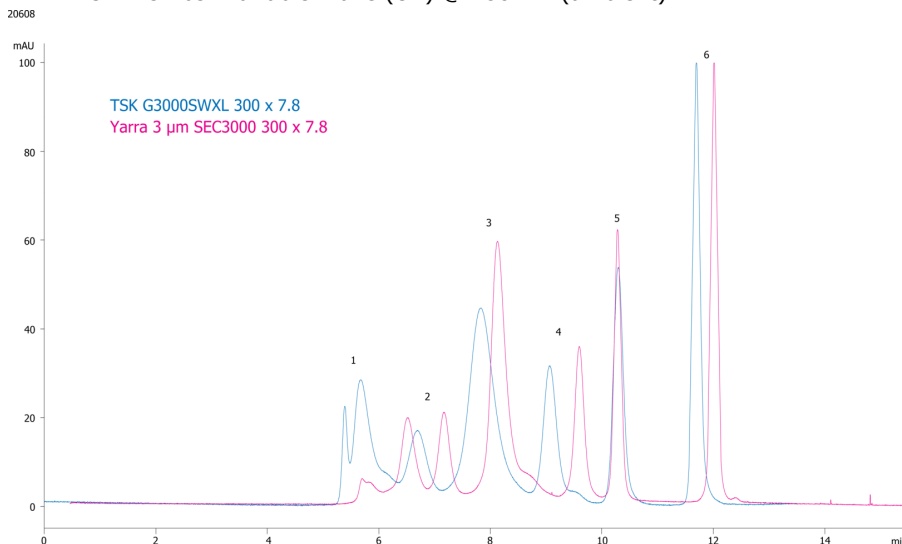

Products used in this application:



ANALYTES:

- 1** Thyroglobulin
Retention Time: 6 min
- 2** IgA
Retention Time: 6.7 min
- 3** IgG
Retention Time: 7.7 min
- 4** Ovalbumin
Retention Time: 9.2 min
- 5** Myoglobin
Retention Time: 9.8 min
- 6** Uridine
Retention Time: 11 min

